

# Concentrated and Ready to Use General Surface Sanitiser



**The General Surface Sanitiser possesses dual action cleaning and disinfecting properties for all hard surfaces. With regular use, the sanitiser offers protection from both bacteria and enveloped viruses.**

It is tested according to BS EN 1276, BS EN 1040 and EN 14476:2013 + A2:2019 Annex A and is proven to have broad range efficacy against bacteria, including *E. coli*, *Paeruginosa*, *E. hirae* and *S. aureus*, and viruses, including all coronaviruses and SARS-CoV-2 (the virus which causes COVID-19).

To use the Concentrated General Surface Sanitiser as an anti-viral, dilute it with water at a ratio of 1:9. As way of example, to produce 750ml of cleaning solution, simply mix 75ml of concentrated sanitiser with 675ml of water. Then spray onto a surface, wait five minutes and wipe clean with a cloth.



CellPath can also provide a ready to use mixture in refillable trigger operated spray bottles.

The General Surface Sanitiser is safe for use on all types of hard surfaces within the laboratory such as benchtops, sinks, tiles, floors, desks and doors/handles.

**Warning:** Always use biocides safely and read the label and product information before use.



## KEY BENEFITS

- 1 Can be used as both an anti-viral and anti-bacterial
- 2 Independent testing confirms efficacy against all enveloped viruses as defined in EN 14476:2013 + A2:2019 Annex A, including all coronaviruses and SARS-CoV-2 (the virus which causes COVID-19)
- 3 Biocidal effectiveness in compliance with the following standards: EN1276, EN1650, EN13697 and EN 14476+A2
- 4 Readily biodegradable with low toxicity, minimising environmental impact
- 5 Contains no quaternary ammonium salts
- 6 Manufactured in accordance with recognised international standards, ISO 9001 and ISO 14001, to ensure consistent quality of product

Item Code	Description	Quantity
XTA-0400-63A	Concentrated General Surface Sanitiser – 5L	2 x 5 Litre
XTA-0401-50B	Ready To Use General Surface Sanitiser - 750ml (Spray Bottle)	6 x 750ml



**CellPath**   
INNOVATION IN CELLULAR PATHOLOGY

CellPath Ltd

80 Mochdre Enterprise Park,  
Newtown, Powys, SY16 4LE, UK

T: +44 (0)1686 611333

E: sales@cellpath.com

cellpath.com